
HDMI Microscope ADSM201

User Manual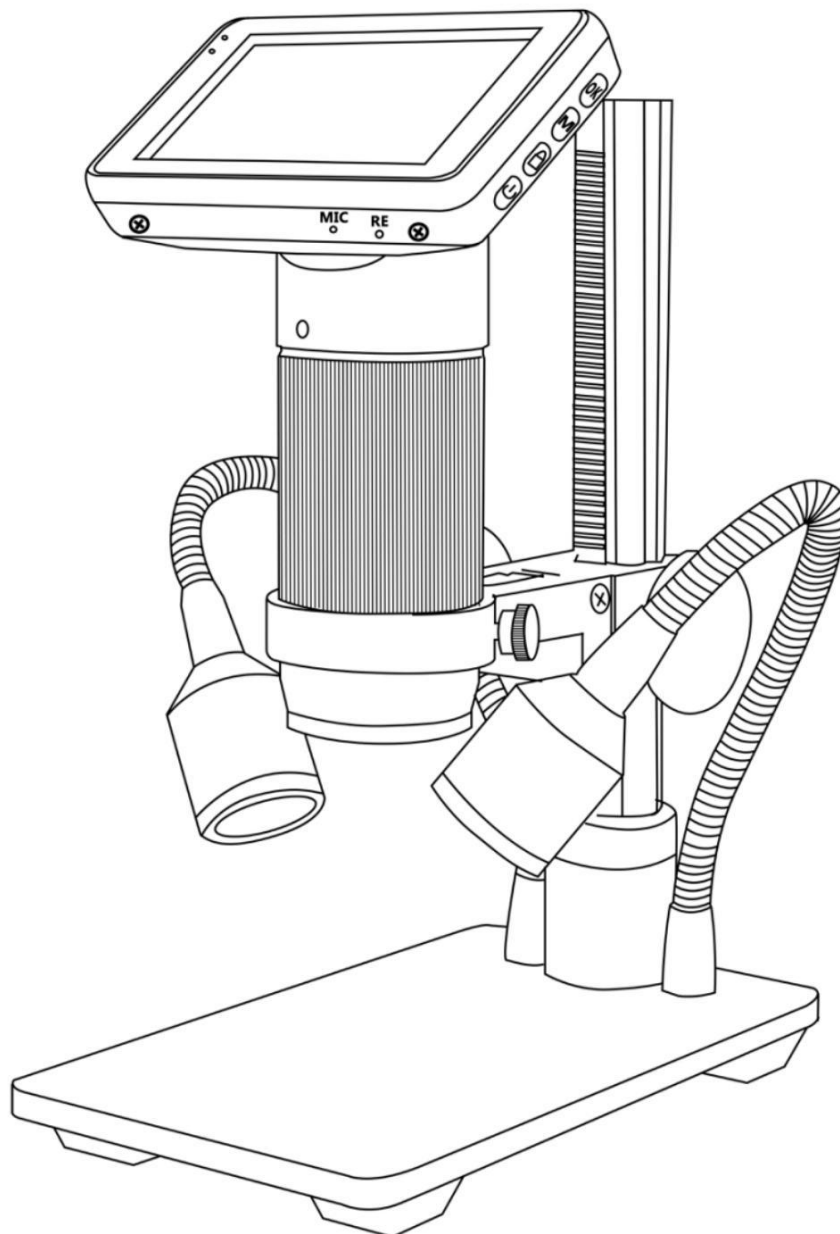

General Rules..... 2

Basic Parameters..... 2

Structure diagram..... 3

Assembly..... 3

General use..... 4

Displaying on an External HDMI Monitor..... 5

USB output mode..... 5

General Rules


ADSM201 is a multifunctional digital microscope. It realized Full HD, high object distance, multifunction, multi-output at the same time. It can easily meet the many different requirements from many industries, such as industrial overhaul and teaching demonstration. And it's appearance design combines many ergonomic consideration, provides users better experience. We sincerely wish this product can release your eyes and improve your work quality. Thanks for your support.

- ADSM201 is safe and low-maintenance, as long as common sense is employed.
- The assembly includes small parts , particularly the thumbscrews used to hold the optical stack in place. Do not allow unsupervised access by infants or toddlers.
- Your ADSM201 is not intended for field use . It is not waterproof, and should not be exposed to extreme temperatures. Treat as you would any other electronic device, such as a laptop or tablet.
- The surface of the lens at the bottom of the ADSM201 should be cleaned whenever you start to notice a degradation in image quality. Use a few drops of lens-cleaning fluid on lens tissues or use a microfiber cleaning cloth.
- The rest of your ADSM201 can be cleaned with a damp cloth.

Basic Parameters

Image sensor	3 Mega Pixels HD Sensor
Output	HDMI&USB
Video format	MOV
Magnification Ratio	Up to 300X(HDMI monitor 22 inch)
Photo resolution	12M
Photo format	JPEG
Focus Range	10mm to 150mm
Frame Rate	Max 30f/s under 600 Lus Brightness
Digital zoom	4X
Storage	Micro-SD card, up to 32G
PC support	Windows XP/7/8
Power source	5V DC
Stand size	14.5cm*7.9cm*16.5cm
Monitor size	3 inch
Accessories	Stand*1, Adapter*1, USB Cable*1, HDMI Cable*1
Packing data	24*18.5*8.3cm/1.2kg

Structure diagram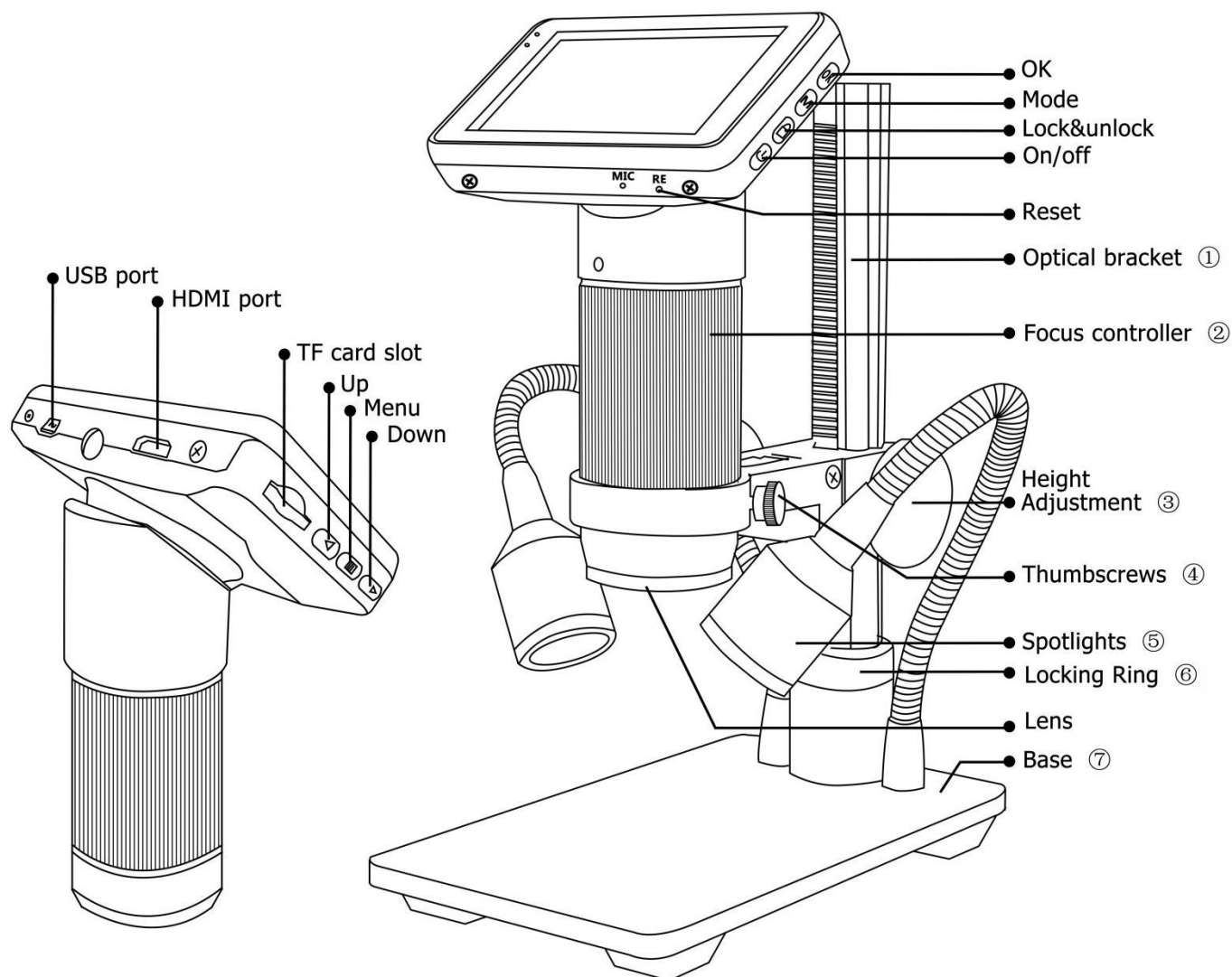

Assembly

1. Bend the spotlights down(5), so that they are out of the way.
2. Screw the locking ring(6) all the way on to the optical bracket(1).
3. Screw the optical bracket(1) on to the base(7). Once you have done so, it is likely the bracket will not be aligned over the base(7). Unscrew it until it is aligned, and then turn the locking ring(6) clockwise until it is tight. (After doing so, it will still be possible to hold the base and turn the bracket if sufficient force is applied, but it will not spin freely.)
4. Loosen the small thumbscrews(4) on the optical bracket(1).
5. Slide the microscope in to the ring of the optical bracket about 3/4" (2 cm), so that the ridged section of the microscope is entirely above the ring.

6. Tighten the thumbscrews(④) to hold the microscope in place. If done correctly, you should be able to turn the focus controller(②) of the microscope freely, while the rest of it remains in place.

7. Remove the protective film from the display at the top of the microscope.

General use

1. Connect the camera to the power supply via USB cable. Also connect spotlights to the power supply.

2. Press the "ON/OFF" button on the monitor. (Sometimes the camera will switch on automatically when it is connected to the power supply)

3. Turn on the spotlights, adjust the "height adjustment" and the "Focus controller" to get the image that you want.

4. Use the "UP/Down" to do digital zooming.

*To set the menu

1. Press "Menu" until it shows the menu on the built-in screen.

2. Press "UP/DOWN" and "OK" choose what to set.

3. Press "Menu" again after you finish setting.

*To switch mode

1. Press "Mode" shortly.

2. Icons :

: standby mode; : capture mode;

 ●: record mode; : image playback mode

: video playback mode (If the files are videos, press "OK" to play the video file)

 ● : Lock the file which is being recording

*To capture pictures.

Press "Mode" shortly to get into the Capture mode. Press "OK" to take pictures.

* To record a video

1.Press "Mode" shortly to get into the Standby mode. Press "OK" to start recording.

2.Press "OK" again to stop recording.

Displaying on an External HDMI monitor


- First, make sure that the monitor connecting to the microscope is literally HDMI monitor. Otherwise it won't perform as good as we expect.
- Second, connect the microscope to the HDMI monitor, also, connect it to the power supply. Because there is no built-in battery in the device, so make sure the microscope is always in charging when you use it.
- Third, turn on the microscope and adjust the distance. Then put the object on the table and adjust the focus controller.
- It won't show anything on its own built-in screen when it is connecting to other monitors. So it's normal that when you use it as HDMI output, the 3 inch screen keeps being dark.
- You still can take pictures and save them if you have a memory card inside the microscope. But you can't record videos in this output mode.

USB output mode

- * The software we provided can only work on Windows XP/7/8 systems.
- * Make sure your computer USB port provided powerful and stable enough electricity. Some laptops may not work well in their condition, please try some desktops USB port, or even the USB port on the back side of the desktop.
- * At this USB output mode, maximum resolution ratio can only get to 1280*720P

Steps:

- 1.Connect the microscope to the computer.
- 2.It will show 2 options, PC camera & Mass storage on the screen of the microscope. Use "Down" button and "OK" button to choose "PC camera" .
- 3.Open the software, it will show the Storage path after the software's name, for example: AMCap-C:\Users\H.chao\Desktop.
- 4.√Select the microscope name in "Devices"
- 5.To take a picture:
 - 1)Aim the storage path of the image file: Capture→Still→Folder→target file.
 - 2)basic settings: Options→Preview√;Capture→Still→Enable√.
 - 3)Take a picture: Capture→Still→Snap.(Or press "Enter" on your keyboard.)
- 6.Record a video:
 - 1)Aim the Storage path of the video file: File→Set capture file→target file
 - 2)Record a video: Capture→start capture→OK→stop capture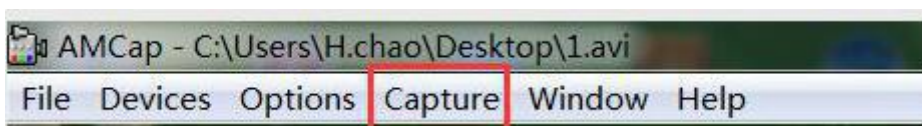

Warranty		
User' s name :	Address :	
Phone number :	Post code :	email :
Model name :	Purchase date :	
Fault feedback :		
Date :	Fault :	
Notice :		
Our warranty doesn' t cover:		
1	If users can not provide the purchase proof or the Warranty.	
2	If users use it in an improper environment, such as incompatible power supply, high temperature environment etc.	
3	If the fault caused by accident, oversight, misoperation or natural disasters etc.	
4	If the fault caused by people who are not belong to authorized organizations of our company during repacking, repair, dismantle, Or if users repack, repair or dismantle the device not following our advice.	
5	Out of warranty period.	

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